

CapTite[™] μCartridge Packing Instructions

CapTite cartridges are easily packed using your packing material and membrane of choice. And when you need to change to a different packing, just disassemble and clean the cartridge and reassemble with your new packing.

- 1. Remove both end caps from cartridge body, remove any old membranes and clean all parts (if applicable).
- 2. Cut or punch two membrane circles ~5mm diameter (~3/16").
- 3. Insert a membrane circle on one end of the cartridge.
- 4. Install an end cap over the membrane and hand tighten.
- 5. Load the open end of the cartridge with a slurry pack to just past the upper membrane level (use syringe or similar to fill).
- 6. Attach a vacuum line to the lower port and evacuate the liquid in the cartridge (ensure the vacuum has a filter to capture the water). Alternatively, a one-piece fitting syringe can also be used for this application.
- 7. Check the level of the packing to ensure it is near but not above the upper membrane level, remove or add additional packing if necessary.
- 8. Insert the second membrane.
- 9. Install an end cap over the top membrane and hand tighten.
- 10. Connect to capillary tubing using the one-piece fittings. Ensure capillary tubing does not extend through the membrane material.

